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
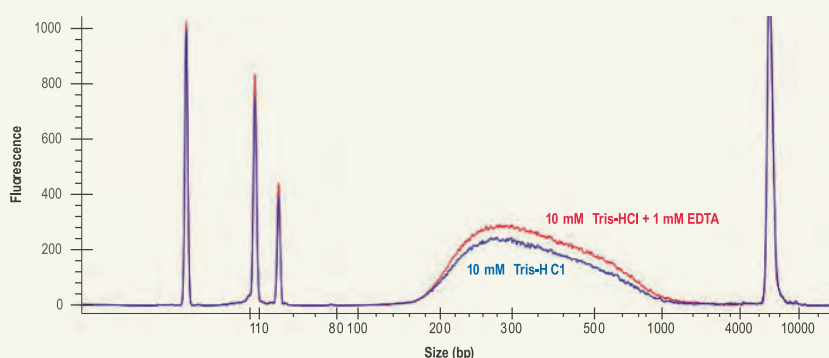
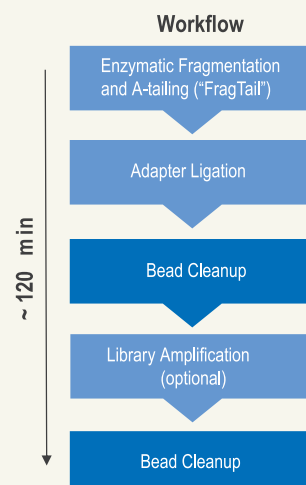
The KAPA EvoPlus chemistry enables reproducible enzymatic fragmentation across DNA diluted in different buffer types.

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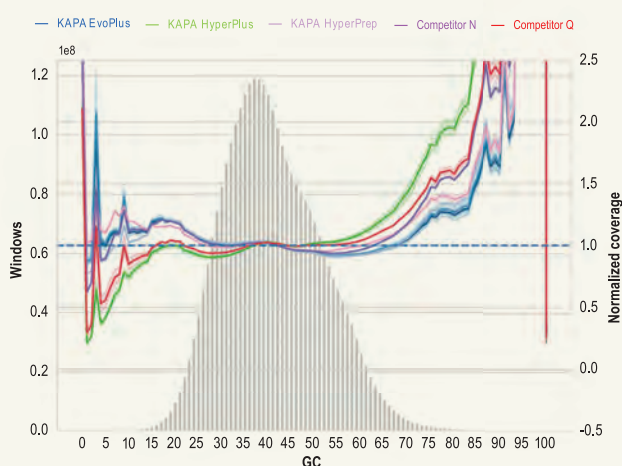
Bias associated with enzymatic fragmentation in the KAPA EvoPlus workflow had no impact on overall coverage depth and uniformity, or GC bias—which were virtually identical for KAPA EvoPlus and KAPA HyperPrep workflows.

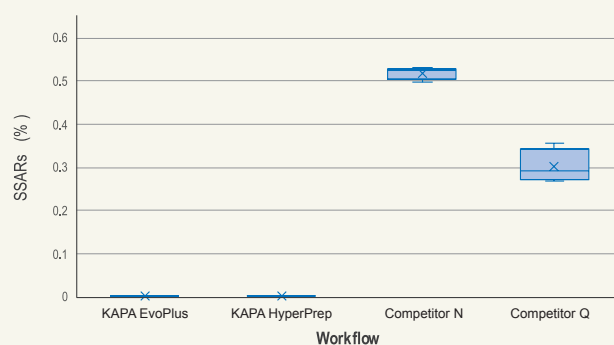優化酵素降低 sequencing artefacts

Improved sequencing performance – reduction of strand-split artefact reads (SSARs).

貨號	品名	包裝
09420037001	KAPA EvoPlus Kit	24 rxns
09420053001		96 rxns
09420339001		384 rxns
09420428001		96 well plate

For research use only, not for clinical purpose.

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

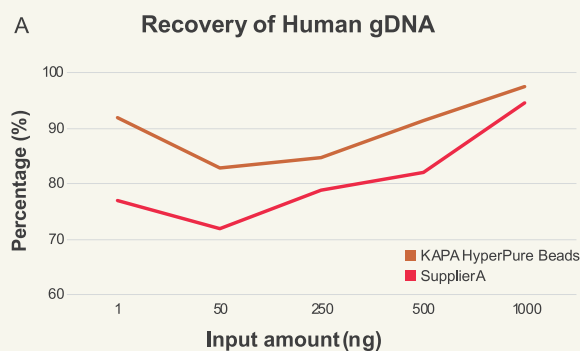
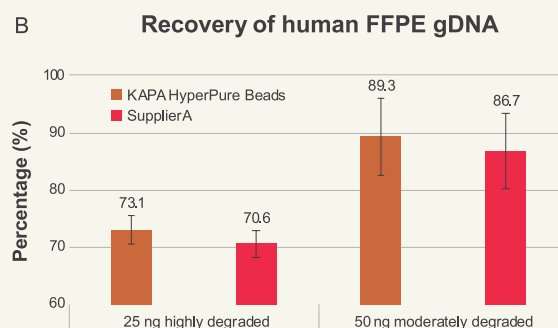
KAPA HyperPure Beads and Supplier A beads were used at a 3X ratio to clean up DNA inputs of various amounts. Recovery was measured using the Qubit Fluorometer 3 dsDNA HS Assay Kit before and after cleanup.

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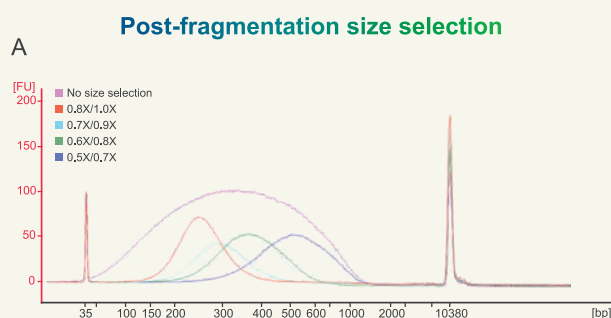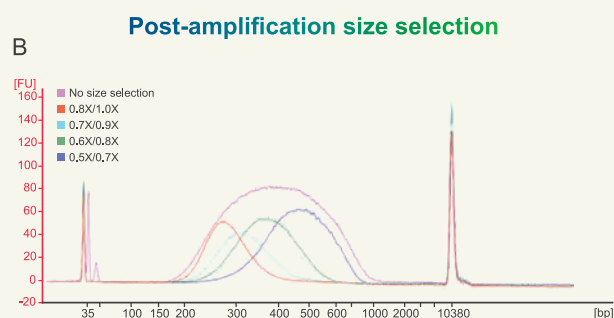High-quality hgDNA was mechanically fragmented with a Covaris E220 Focused Ultrasonicator using conditions optimized to yield a mode fragment length of 250 – 400 bp.

Libraries were PCR amplified and final amplified libraries were subjected to size selection.

貨號	品名	包裝
08963835001	KAPA HyperPure Beads	5 mL
08963843001		30 mL
08963851001		60 mL
08963878001		4 x 60 mL

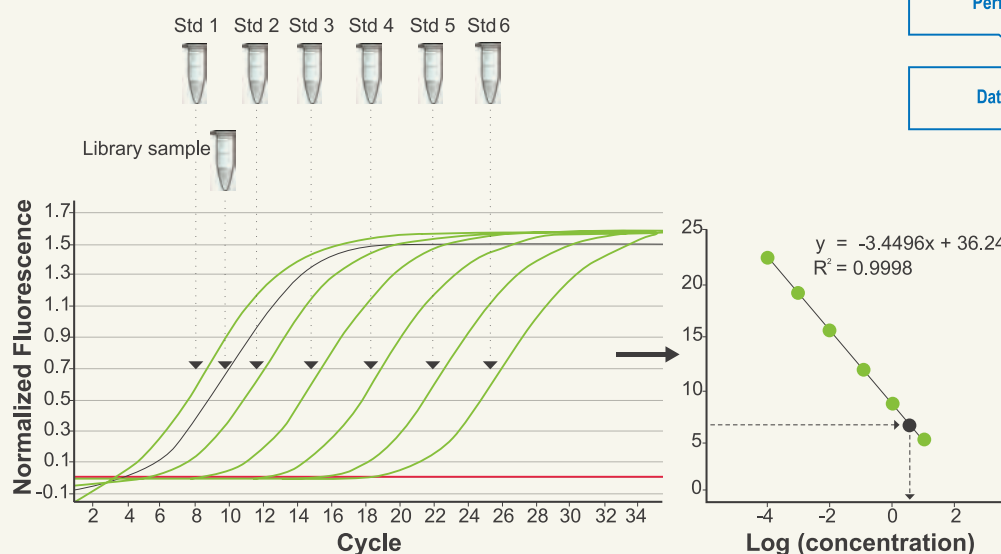
For research use only, not for clinical purpose.

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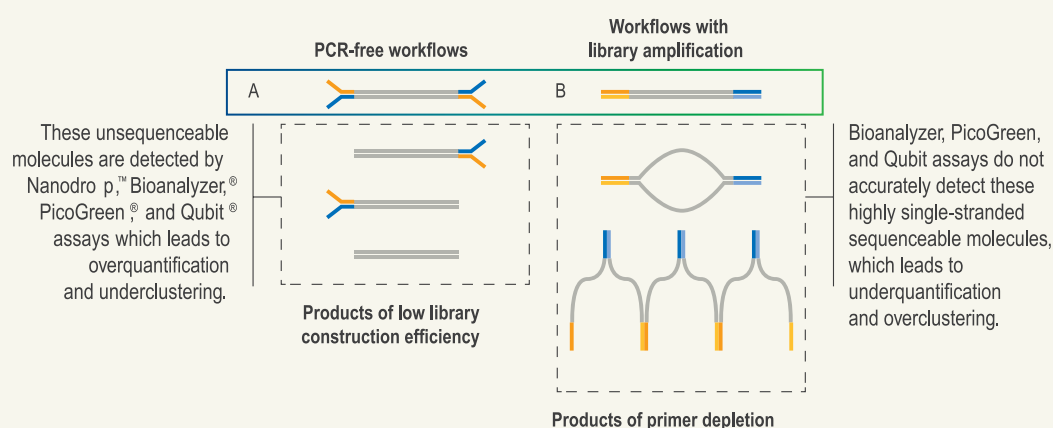
for research use only, not for clinical purpose

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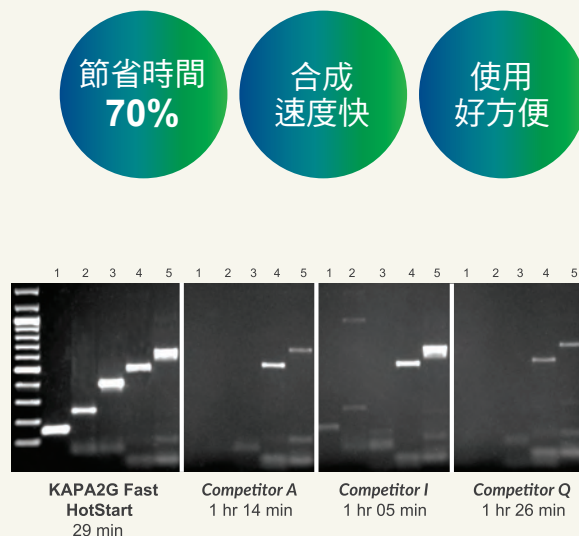
Reactions (25 μ L) contained 5 ng human genomic DNA and 0.5 units (KAPA2G Fast HotStart and Competitor I) or 0.625 units (Competitor A and Q) enzyme. The extension (72°C) was 1 sec/cycle for KAPA2G Fast HotStart and 60 sec/cycle for competitor enzymes.

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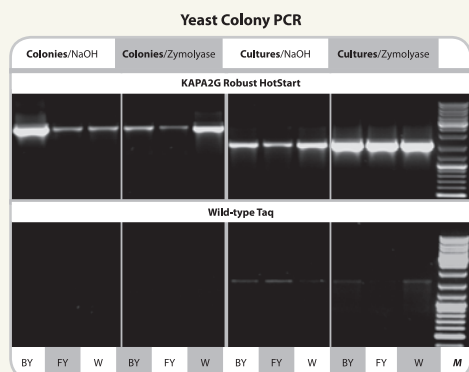
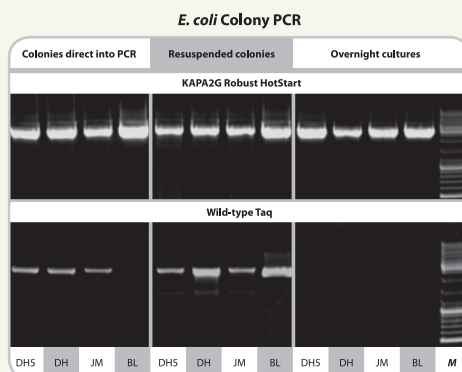
Amplification of a 2.5 kb (left) or 1.6 kb (right) fragment of the GSH1 gene from three commonly used *S. cerevisiae* strains (BY4742, FY23 and W303). A standard 3-step cycling profile (35 cycles), with 30 sec/kb extension time was used for KAPA2G Robust HotStart reactions. For wild-type Taq, the extension time was 1 min/kb.


Amplification of a 2.7 kb insert from four commonly used *E. coli* strains (DH5a, DH10B, JM109 or BL21). A standard 3-step cycling profile (35 cycles), with 30 sec/kb extension time was used for KAPA2G Robust HotStart reactions. For wild-type Taq, the extension time was 1 min/kb.

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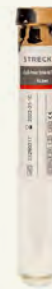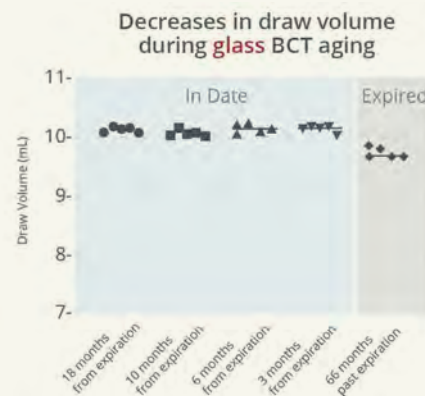
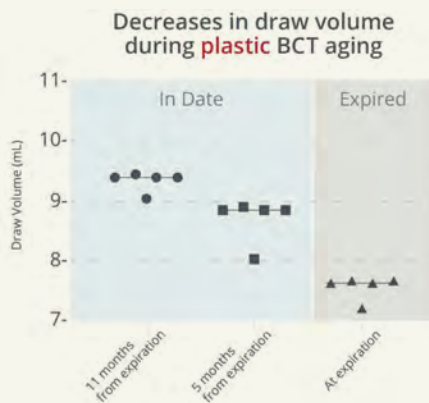
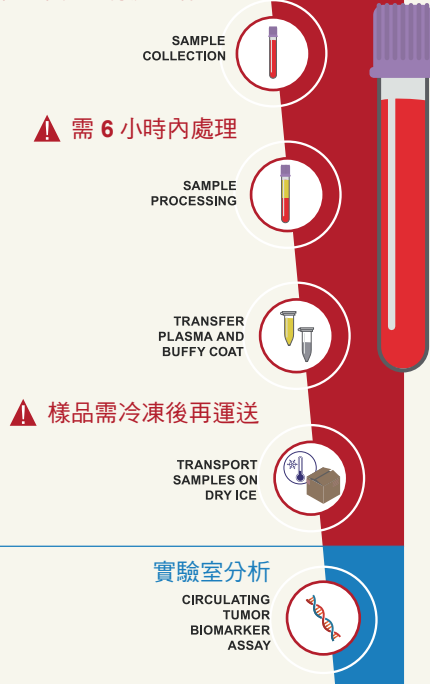
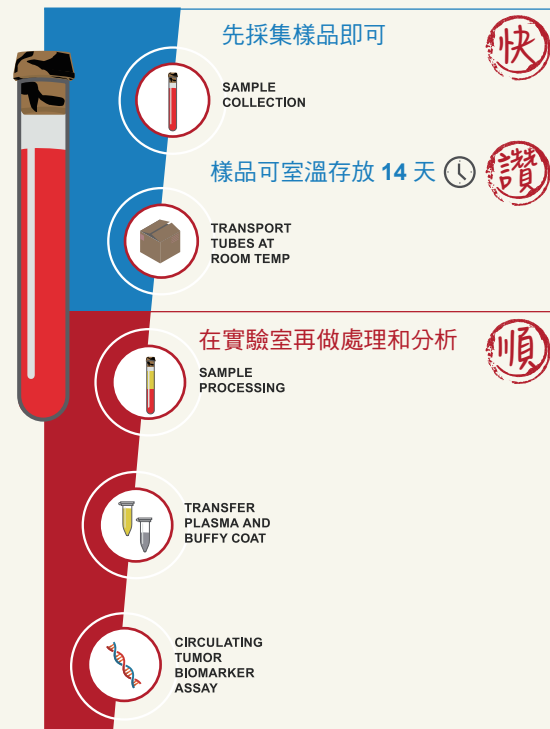
RNA extraction Selection guide
Acid-guanidium-phenol based reagents
例如：TRIzol®, TriPure, RNAzol, TRI Reagent


貨號	品名	包裝	促銷
R1051	Quick-RNA™ MicroPrep Kit	200 Preps	買大送小 R1050
R1055	Quick-RNA™ MiniPrep Kit	200 Preps	買大送小 R1054
R1058	Quick-RNA™ MiniPrep Plus Kit	200 Preps	買大送小 R1057
R2052	Direct-zol™ RNA MiniPrep	200 Preps	買大送小 R2050
R2053	Direct-zol™ RNA MiniPrep 含 200 ml TRI reagent	200 Preps	買大送小 R2051
R2062	Direct-zol™ RNA MicroPrep	200 Preps	買大送小 R2060
R2063	Direct-zol™ RNA MicroPrep 含 200 ml TRI reagent	200 Preps	買大送小 R2061
R2072	Direct-zol RNA Miniprep Plus	200 Preps	買大送小 R2070
R2073	Direct-zol RNA Miniprep Plus 含 200 ml TRI reagent	200 Preps	買大送小 R2071

Bisulfite Conversion-The Gold Standard

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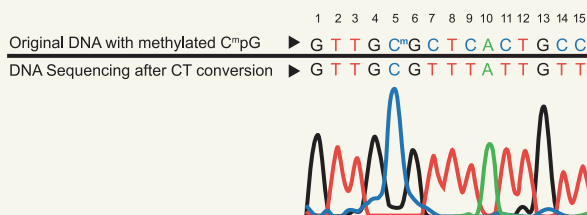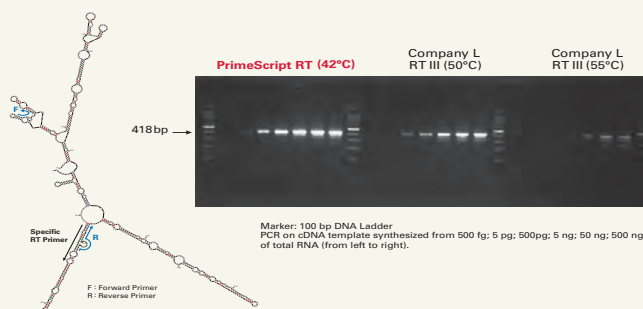
Mir-X™ miRNA First Strand Synthesis Kit

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

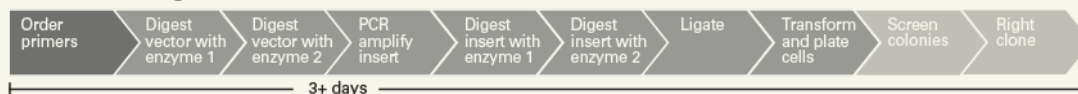
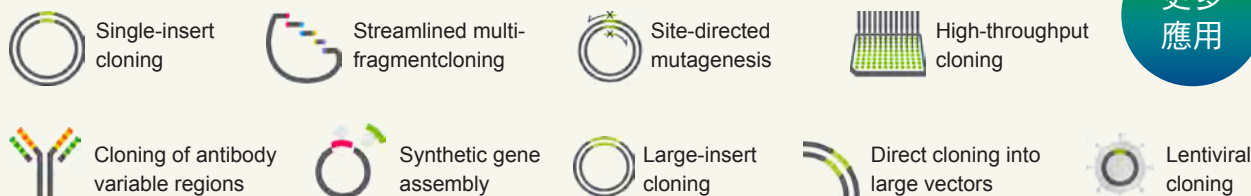
CLONTECH Retro-X / Lenti-X Concentrator

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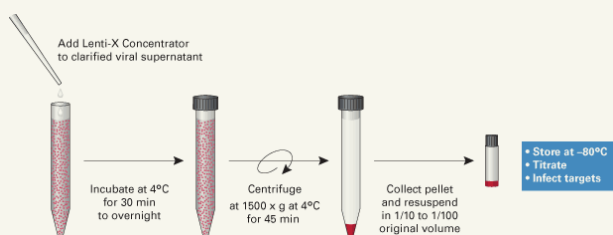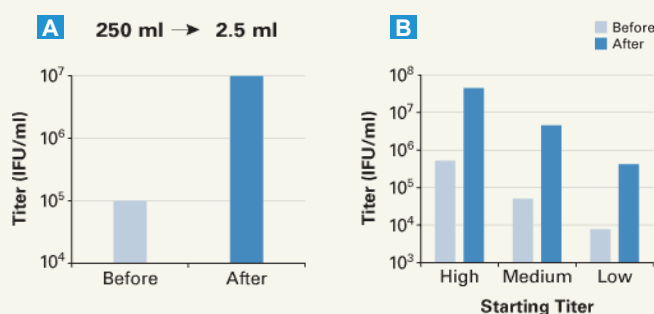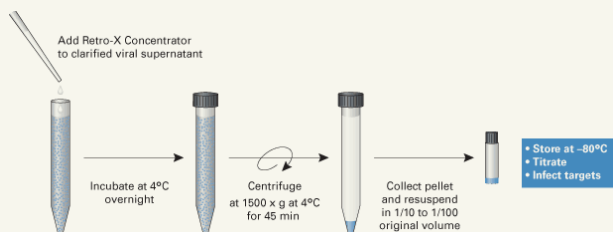
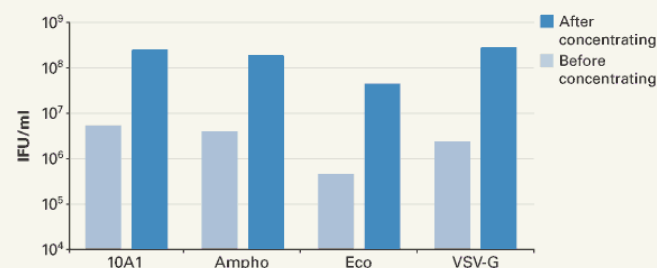
References

- Morelli, Kathryn H et al. "An RNA-targeting CRISPR-Cas13d system alleviates disease-related phenotypes in Huntington's disease models." *Nature neuroscience* vol. 26,1 (2023): 27-38. doi:10.1038/s41593-022-01207-1
- Kim, Hark Kyun et al. "Engineered small extracellular vesicles displaying ACE2 variants on the surface protect against SARS-CoV-2 infection." *Journal of extracellular vesicles* vol. 11,1 (2022): e12179. doi:10.1002/jev2.12179
- Sun, Tianliang et al. "ZNR3 and RNF43 cooperate to safeguard metabolic liver zonation and hepatocyte proliferation." *Cell stem cell* vol. 28,10 (2021): 1822-1837.e10. doi:10.1016/j.stem.2021.05.013
- Alam, Sk Kayum et al. "ASCL1-regulated DARPP-32 and t-DARPP stimulate small cell lung cancer growth and neuroendocrine tumour cell proliferation." *British journal of cancer* vol. 123,5 (2020): 819-832. doi:10.1038/s41416-020-0923-6
- Abdeladhim, Maha et al. "Engineered ovalbumin-expressing regulatory T cells protect against anaphylaxis in ovalbumin-sensitized mice." *Clinical immunology (Orlando, Fla.)* vol. 207 (2019): 49-54. doi:10.1016/j.clim.2019.07.009

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TFF-Easy Tangential Flow Filter for EV concentration

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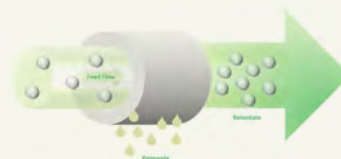
貨號	產品	包裝	Pore size
HBM-TFF/1 HBM-TFF/5	TFF-Easy - tangential flow filtration EV concentrator	1 concentrator 5 concentrators	5 nm
HBM-TFF-EVs-S HBM-TFF-EVs-L	TFF-EV- tangential flow filter for EV purification	1 concentrator	50 nm
HBM-TFF-MV	TFF-MV- tangential flow filter for large EV separation and recovery	1 concentrator	200 nm

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Virus concentration reagent- PEG-it™

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Endothelial Cell Growth Supplement (ECGS)

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
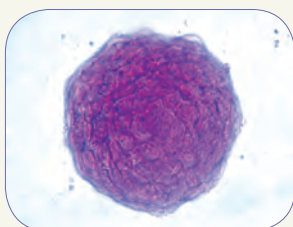
Alkaline Phosphatase staining of Mouse ES cells.

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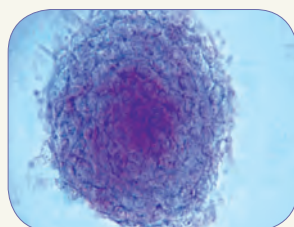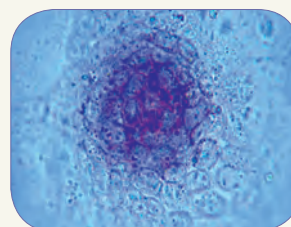
Compartment Protein Extraction Kit

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貨號	品名	包裝
2145	Compartment Protein Extraction Kit	5 grams tissues or 125 million cells

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ProteoExtract®

Native Membrane Protein Extraction Kit

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
BugBuster® Protein Extraction Reagent

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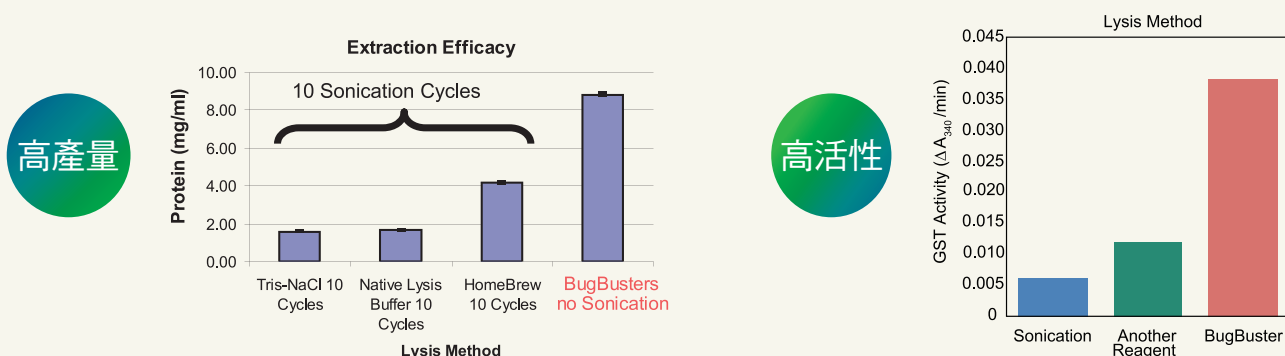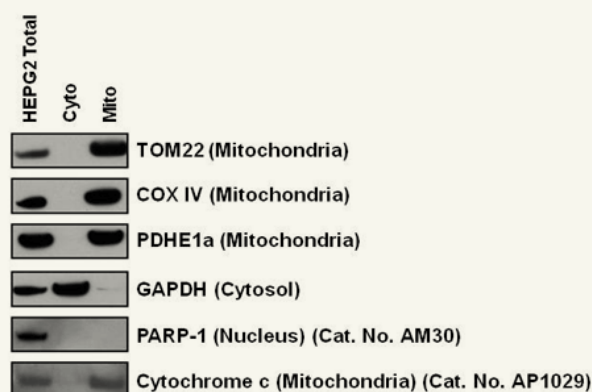
Protein Extraction Reagents Application Guide

Cat No. cell type product			Starting Material		Applications							
			Total Culture	Cell Pellet	HT Compatible	Analysis						Purification
						1D PAGE	2D PAGE	IEF	MS	WB	Activity Assay	
E-coli												
71456	BugBuster Master Mix			●	●	●	●		●	●	●	
70584	BugBuster Protein Extraction Reagent			●		●	●		●	●	●	
70922	BugBuster HT Protein Extraction Reagent			●	●	●	●		●	●	●	
70923	BugBuster (Primary amine-free) Extraction Reagent			●		●	●		●	●	●	
70921	BugBuster 10X Protein Extraction Reagent			●	●	●	●		●	●	●	
71092	PopCulture Reagent		●		●	●			●	●	●	
Yeast												
71186	YeastBuster Protein Extraction Reagent			●		●			●	●	●	
Insect												
71009	CytoBuster Protein Extraction Reagent			● ⁺		●	● [*]	● [*]		●	●	
71187	Insect Protein Extraction Reagent		●		●	●			●	●	●	
Lysis and Extraction Enhancement												
71110	Gram-negative bacteria (E-coli)	rLysozyme Solution	●	●	●	●			●	●	●	
71230		rLysozyme Bioprocessing Reagent	●	●	●	●			●	●	●	
71412	Gram-positive bacteria (E-coli)	Chicken Egg White Lysozyme Solution	●	●	●	●			●	●	●	
70746	All cells	Benzonase Nuclease	●	●	●	●			●	●	●	

Key : 1D PAGE = One-dimensional Polyacrylamide Gel Electrophoresis
 2D PAGE = Two-dimensional Polyacrylamide Gel Electrophoresis
 * = Salt must be removed before IEF

MS = Mass Spectrometry R = Reporter Assay +=Monolayer
 IEF = Isoelectric Focusing G = Gel Shift

Mammalian Cell Protein Extraction Reagents Application Guide

分類	貨號	品名	包裝
核蛋白	2900	Nuclear Extraction Kit Mammalian cells	100 Rxns
膜蛋白	444810-1KITCN	ProteoExtract® Native Membrane Protein Extraction Kit	20 Extractions
	71772-3CN	ProteoExtract® Transmembrane Protein Extraction Kit	20 Extractions
粒線體	QIA88-1KITCN	ProteoExtract® Cytosol/Mitochondria Fractionation Kit	100 Extractions
Subcellular Protein	539790-1KITCN	ProteoExtract® Subcellular Proteome Extraction Kit	20 Extractions
	2145	Compartment Protein Extraction Kit	Kit (5g tissue / 1.25 x 10 ⁹ cells)
Total protein	2140	Total Protein Extraction Kit	Kit (5g tissue / cells)
	71009-4CN	CytoBuster™ Protein Extraction Reagent	250 mL
	20-188	10X RIPA Lysis Buffer	100 mL

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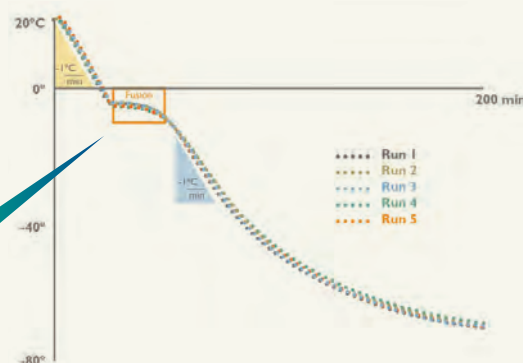
CoolCell LX Cell Freezing Containers 12 孔

貨號	Description	Color	Dimensions
CORNING 432001	CoolCell LX, Cell Freezing Container for 12 x 1ml to 2ml cryogenic vials	Blue	Diam 11.7 X 10.9 cm
CORNING 432002	CoolCell LX, Cell Freezing Container for 12 x 1ml to 2ml cryogenic vials	Green	Diam 11.7 X 10.9 cm
CORNING 432003	CoolCell LX, Cell Freezing Container for 12 x 1ml to 2ml cryogenic vials	Orange	Diam 11.7 X 10.9 cm
CORNING 432004	CoolCell LX, Cell Freezing Container for 12 x 1ml to 2ml cryogenic vials	Pink	Diam 11.7 X 10.9 cm

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CoolCell FTS30 High-Capacity Cell Freezing Containers 30 孔

貨號	Description	Color	Dimensions
CORNING 432006	CoolCell FTS30, Cell Freezing Container for 12 x 1ml to 2ml cryogenic vials	Blue	Diam 16.5 X 11.5 cm
CORNING 432007	CoolCell FTS30, Cell Freezing Container for 12 x 1ml to 2ml cryogenic vials	Orange	Diam 16.5 X 11.5 cm
CORNING 432008	CoolCell FTS30, Cell Freezing Container for 12 x 1ml to 2ml cryogenic vials	Green	Diam 16.5 X 11.5 cm
CORNING 432009	CoolCell FTS30, Cell Freezing Container for 12 x 1ml to 2ml cryogenic vials	Pink	Diam 16.5 X 11.5 cm

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CoolCell LX Cell Freezing Containers

貨號	Description	Color	Dimensions
CORNING 432005	CoolCell 5ml LX, Cell Freezing Container for 12 x 3.5ml to 5ml cryogenic vials	Blue	Diam 9.5 X 14.5 cm

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