# Liquid Biopsy

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
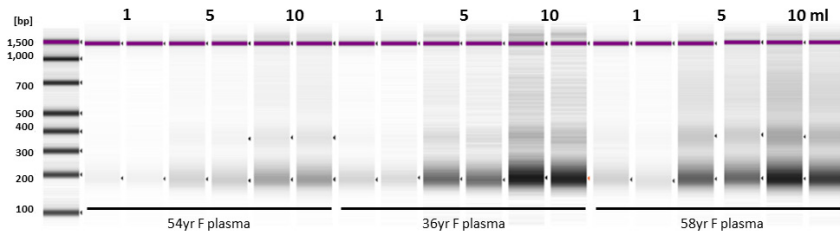
Total DNA is efficiently purified from cell-free biological fluids with the Quick-cfDNA™ Serum & Plasma Kit. Total DNA, including both high and low molecular weight species, purified (duplicates) from human maternal plasma, amniotic fluid and cerebrospinal fluid was analyzed.

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
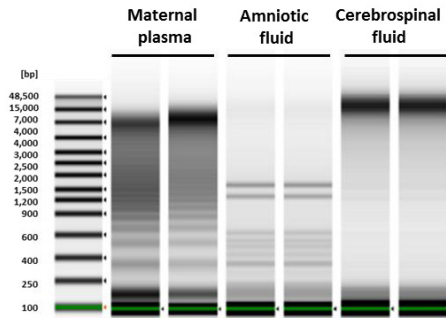
Cell-free DNA recovery scales proportionally with sample input using the Quick-cfDNA™ Serum & Plasma Kit gel image show the linear recovery of cfDNA from human plasma and serum (healthy female donors), as measured.

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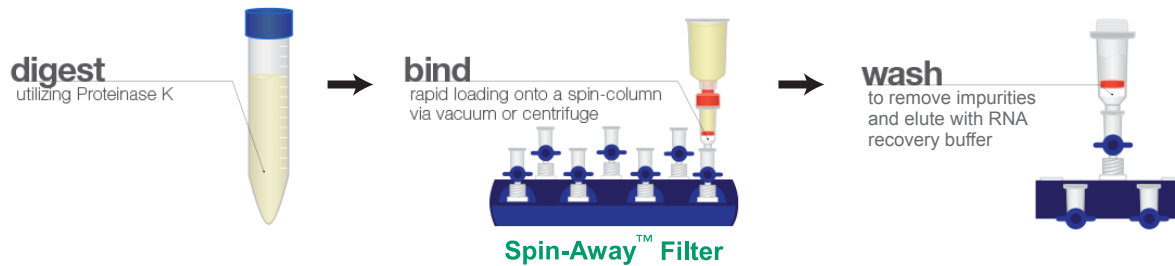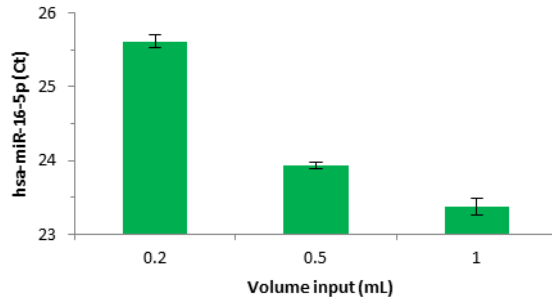
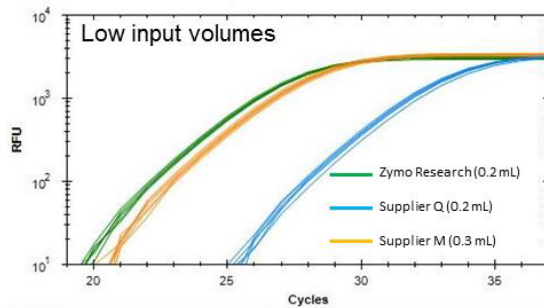
Proportional scale up relative to input volume (from 0.2 ml to 1 ml; 55y male plasma). Triplicates of purified samples used to generate RT-qPCR (BioLine; EPIK miRNA Select Kit, hsa-miR-16-5p assay).

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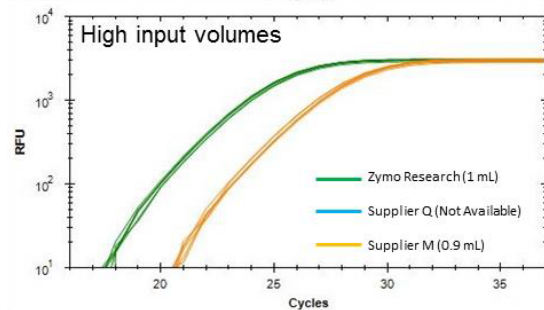
2x and 8x more yields achieved from low and high input volumes, respectively, compared to the next best product from other supplier. Triplicates of purified samples from different kits used to generate RT-qPCR data .

# 血漿游離 DNA (cfDNA) & 外吐小體 (Exosomes)
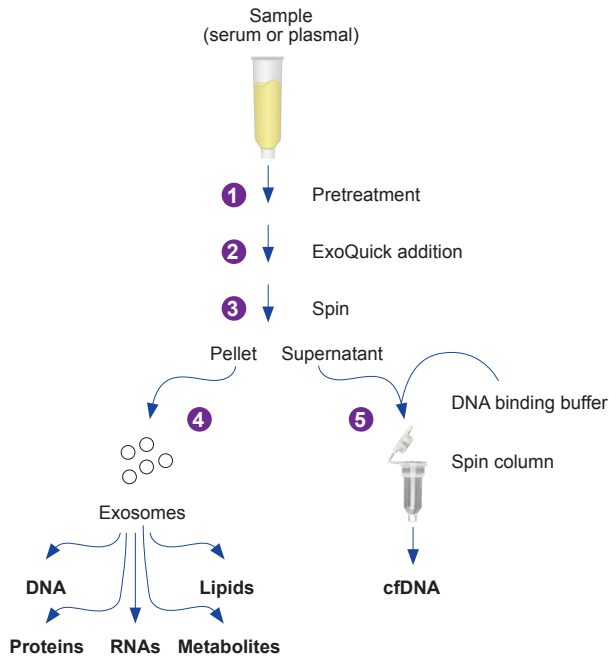
## XCF™ COMPLETE Exosome & cfDNA Isolation Kits

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

### non-exosomal protein 殘留率低

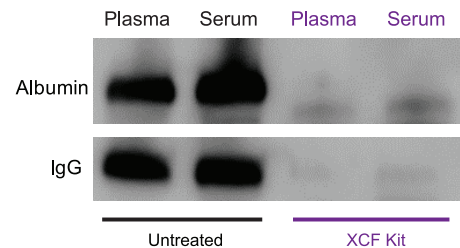The XCF COMPLETE Kit shows low protein carryover. Western blots of co-purifying human albumin (~67kD) and heavy-chain IgG (~50kD) comparing untreated serum and plasma versus the same samples processed using the XCF Kit .

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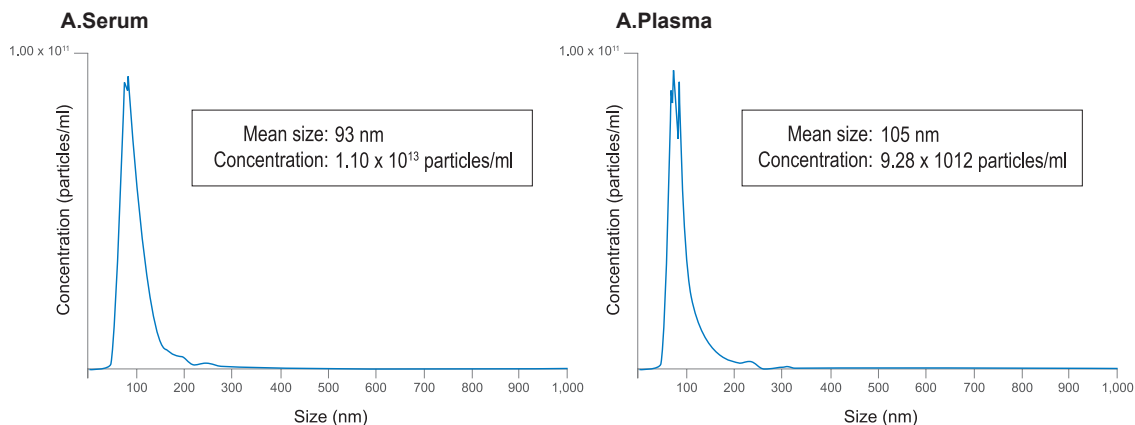
Exosomes prepared using the XCF COMPLETE Kit deliver the expected particle size profile. NanoSight NTA data for extracellular vesicles isolated from 500  $\mu$ l of serum (Panel A) and plasma (Panel B).

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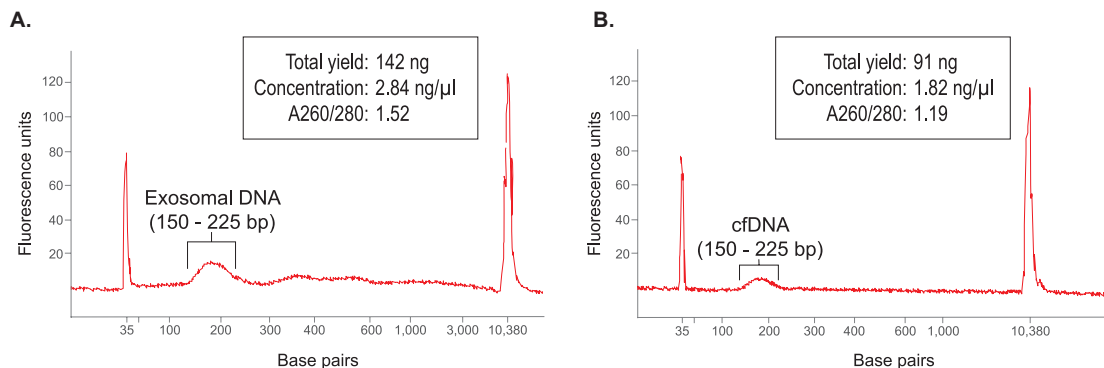搭配 **XCF™ Exosomal DNA Isolation Kit** 也可以萃取出 **exosomal DNA**

cfDNA and exosomal DNA isolated using the XCF COMPLETE Kit are of good quality and yield with the expected size. Agilent Bioanalyzer profiles of the exosomal DNA portion (Panel A) and cfDNA portion (Panel B) of 500  $\mu$ l of serum processed with the XCF COMPLETE Kit. A notable peak around ~166 bp is seen in both cfDNA and exosomal DNA.

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Nucleotide cut-off guidelines for single stranded DNA

Single-Stranded Nucleotide Cut-off (bp)	NMWL	
	3K (40 min)	10K (20 min)
10	98	97
15	96	95
20	94	94
25	96	96

Nucleotide cut-off guidelines for double stranded DNA

Double-Stranded Nucleotide Cut-off (bp)	NMWL		
	30K	50K	100K
137	95	95	—
301	—	94	—
657	—	99	—
1129	—	100	97

Based on >90% retention. Spinning 50 and 100K devices at higher than recommended g-force may result in lower recovery of DNA.

Recommended g-force and spin time for Amicon Ultra devices

NMWL	Maximum G-Force Rating	Spin Time (min) at 25 °C
3K	7,500	40
10K	7,500	20
30K	5,000	10
50K	2,000	20
100K	2,000	10

	Cat. No.	name	size
	純化	血漿游離 DNA (Cell-free DNA; cfDNA) & 血漿游離 RNA (Cell-free RNA; cfRNA)	
 ZYMO RESEARCH <i>The Beauty of Science is to Make Things Simple</i>	D4076	Quick-cfDNA™ Serum & Plasma Kit	50 preps
	R1059	Quick-cfRNA™ Serum & Plasma	50 preps
 SBI System Biosciences	XCF100A-1	XCF COMPLETE Exosome and cfDNA Isolation Kit	20 rxns
	XCF200A-1	XCF Exosomal DNA Isolation Kit	20 rxns
	濃縮	循環腫瘤 DNA ( circulating tumor DNA; ctDNA )	
 MERCK	UFC500324	Amicon Ultra-0.5 Centrifugal Filter Unit ; 3 kDa	24/pk
	UFC501024	Amicon Ultra-0.5 Centrifugal Filter Unit ; 10 kDa	24/pk
	UFC503024	Amicon Ultra-0.5 Centrifugal Filter Unit ; 30 kDa	24/pk
	UFC505024	Amicon Ultra-0.5 Centrifugal Filter Unit ; 50 kDa	24/pk
	UFC510024	Amicon Ultra-0.5 Centrifugal Filter Unit ; 100 kDa	24/pk